



## Integration with Liquid Handling Robots

### A Integration with PerkinElmer Multiprobe



The pictures above show one possibility to integrate the **chemagic Magnetic Separation Module I** (chemagic MSM I) with a PerkinElmer Multiprobe Liquid Handling instrument. The right segment of the Multiprobe is substituted by the horizontal tracking unit of the chemagic MSM I. The gripper is able to place and remove 12 tube racks (Falcon format) or 96 well plates on the tracking unit enabling a fully automated process for high throughput sample preparation.

### B Integration with Tecan Genesis (EVO)



An integration with Tecan instruments works in a comparable manner. The main difference is the position of the chemagic MSM I which in this case is situated under the Tecan platform. A large size RoMa (Robotic Manipulator) of the Liquid Handler enables the positioning of racks or plates on the tracking unit.

*Further Questions?*

**Phone** +49 (0) 2401 805-501

**chemagen** technical  
support

**Mail** [support@chemagen.de](mailto:support@chemagen.de)